© American Scientific, LLC AmSci Stream Table Kit Item # 6600-00

## KIT SET-UP

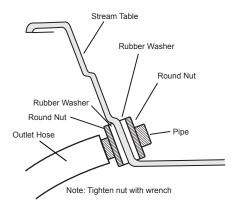


Figure 1

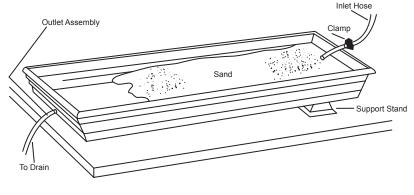


Figure 2

- d. Direct the outlet hose into a sink, drain, or pail. See Figure 2.
- 3. Prepare the inlet hose:
  - a. Attach the second plastic hose to a water faucet, recirculation pump or siphon system.
  - b. Direct the other end of the hose through the adjustable clamp and into the stream table opposite the outlet hose.
  - c. Adjust the flow of water using the adjustable clamp.
- 4. Place the plastic supports under the inlet hose end of the Stream Table so that it is raised higher than the outlet end. See Figure 2.